User Manual 用户手册

EN 耐火安全储存柜
Safety Storage Cabinets

SYSBEL® **SYSBEL IS SAFETY**
安全·源自西斯贝尔

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Thank you for using our products. As a company of caring employee safety and environmental protection, SYSBEL provides high quality products to ensure the safety of your working environment and personnel.

Description

1. The cabinet is made of high-strength steel plates, the surface of the cabinet is Electrostatic powder spraying which makes the surface easy to clean, scratch-resistant, shock resistant, and durable. The inner of the cabinet consists of multiple layers of fireproof material in order to prevent the fire for 90 minutes.
2. When using the safety cabinet in a normal situation, the door of the cabinet can be opened or closed manually; if there is a fire accident, when the temperature reaches 122°F (50°C), the built-in fuse will be connected to the system and close the opened cabinet door automatically in order to protect the item inside the cabinet. There is a damper in the cabinet door which can adjust the force of open and close the door.
3. The Vents comply with EN14470-1. The Vent baffle will automatically close at 158°F (70 ° C) in a fire accident. The ventilation system can be externally connected with an air exhausting pipe or a circulating air filter;
4. When there is a fire accident, the special heat sensitive material in the seals and all the connection areas can expand automatically and fill the seals to keep the inner of the cabinet in a sealed status.
5. The standard cabinet is equipped with 3 adjustable steel shelves, the adjustable distance between each shelf is 32mm, the bearing weight of shelf is 75 kg. The edge of the shelf is 30mm higher than the inner in order to prevent the small vessels from dropping by accident.
6. Each door has a separate cylinder lock to effectively prevent unauthorized use. Reflective warning labels in multiple languages are displayed on the door to stand out under the light of flashlight in the case of fire or power outage. The label complies with ISO 3864 international standard.
7. The detachable leakage-absorbed sump in the bottom is used in conjunction with perforated inserts to effectively absorb the overflow liquid on the shelf. Containers placed on perforated inserts can avoid being soaked in the bottom of the drain sump;
8. The bottom of the cabinet body is equipped with adjustable level support feet, which can keep the balance of the cabinet even on the uneven surface;
9. The cabinet is equipped with an electrostatic grounding conducting port for easy connection of anti-static device
10. The safety cabinet complies with EN 14470-1 and regulation of NFPA30 and OSHA as well. These regulations ensure the properties and structure requirements of the cabinet, providing safe storage space for flammable liquids to reduce the fire risks.

TÜV Certification

SE860230

TÜV Certification

SE830450

TÜV Certification

SE890450

TÜV Certification

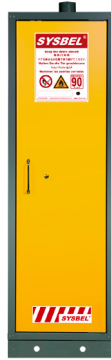
TÜV The actual test passed the sample

Comply with EN, NFPA30 and OSHA standards

Under Counter Cabinet

SE890230

SE490300

WAB005

Safety Storage Cabinets												
Fire rating (minutes)	Door type	Color	Standard	Ext dimension (H×W×D/cm)	Packing dimension (H×W×D/cm)	Adjustable Shelves (Pcs)	Shelf Loading Weight (Kgs.)	Capacity of bottom tank (L)	Diameter of air outlet connection (mm)	Can store a single largest container (L)	Volume (gallons /liter)	Model
30	Double Door	Yellow	EN14470-1	204×116×62.5	218×123×74	3	75	37	100	33	45/170	SE830450
60	Single Door			204×60×62.5	218×67×74		75	18.5		16.5	23/87	SE860230
90	Single Door			204×60×62.5	218×67×74		75	18.5		16.5	23/87	SE890230
90	Double Door			204×90×62.5	218×97×74		75	30		27	30/114	SE490300
90	Double Door			204×116×62.5	218×123×74		75	37		33	45/170	SE890450

Safety Under Counter Cabinet													
Fire rating (minutes)	Door type	Color	Standard	Ext dimension (H×W×D/cm)	Packing dimension (H×W×D/cm)	Adjustable Shelves (Pcs)	Shelf Loading Weight (Kgs.)	Capacity of bottom tank (L)	Diameter of air outlet connection (mm)	Can store a single largest container (L)	Volume (gallons /liter)	Model	
90	Single Door	Yellow	EN14470-1	61.5×56×52	75.5×63×63.5	1	75	11	75	9.5	10/38	SE490100	
90	Single Door			75×60×52	89×67×63.5		75	12.5		75	11	12/45	SE490120
90	Double Door			61.5×86×52	75.5×93×63.5		75	19		75	17	17/64	SE490170
90	Double Door			63×110×59	77×117×70.5		75	32		75	28.5	19/72	SE490190

Moving wheel					
Product Name	Loading Weight(Kgs.)	Adaptation cabinet type		Model	Details
Moving wheel	600kg	SE490300. SE830450. SE890450		WAB005 (One group)	Convenient for the cabinet to move indoors

*Selection tips: The colors of this series of products are yellow, blue, white, etc., only the color of the cabinet door is different, the yellow series SKU is as shown in the picture above, the blue series please add "-B" after the SKU in the above table, and the white series For series, please add "-W" after the SKU in the above table

i 使用说明 (Instructions for use)

1. 按照图 1 确认产品外形尺寸 (mm) 和功能点位, 选用合适的工具来搬运到适合放置的区域; 1.
Ensure the product overall dimension(mm) and functional spots according to picture 1, choose the proper tool to help remove the cabinet to the proper area.

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Ensure the level of the cabinet after placing the cabinet, or the incline of the cabinet will lead the door of the cabinet to open and close difficultly. If the cabinet is not level, you can use the No.1 and No.2 of picture 2 to adjust the cabinet. The adjust method is using a professional lifting tool to lift the cabinet 50mm above the ground, and fix the No.1(4pcs) onto the No.5 in a vertical direction, lay the No.2 under the opposite ground of No.5, and adjust No.1 to reach the level of the cabinet. After the cabinet is placed in the level, install the No.4 on the cabinet body through No.3.

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The opening and closing of the lock on the cabinet door: the lock and handle are marked in picture 3. When opening the door, you need to insert the key and rotate clockwise two circles, pull the handle by hand at the same time. When closing the door, you need to turn the key anticlockwise and pull out the key.

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The adjustment of shelf: The standard cabinet is equipped with 3 adjustable steel shelves, the adjustable distance between each shelf is 32mm. Users can adjust the height of each shelf according to the actual need, the adjustment method: as the picture 4 shows, you should remove all the items on No.3 firstly, and then move the No.3 out of the cabinet, use the screwdriver to dismantle two No.1 (8 tapping screw) on the left and right of the No.2, after adjusting the layer to the ideal height, you can use a cross point screwdriver to fix No.2 through No.1 (8 screws) on the left and right side boards, the screws on the left and right side need to be tightened, then place the No.3 on No.2.

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The clean of the leakage liquid in the leakage-absorbed sump: If there is a small amount of leaked liquid on the adjustable shelf, please remove all the stored items on shelf, using the corresponding adsorbents for adjustable shelf clean. If there are any leaked liquids in the bottom leakage-absorbed sump, please try to clean out according to the picture 5, empty items stored on the No.1 firstly, remove the No.1 out of the cabinet, and then use the corresponding adsorbents to clean No.2. After completing the cleaning, follow the steps to put in No.1.

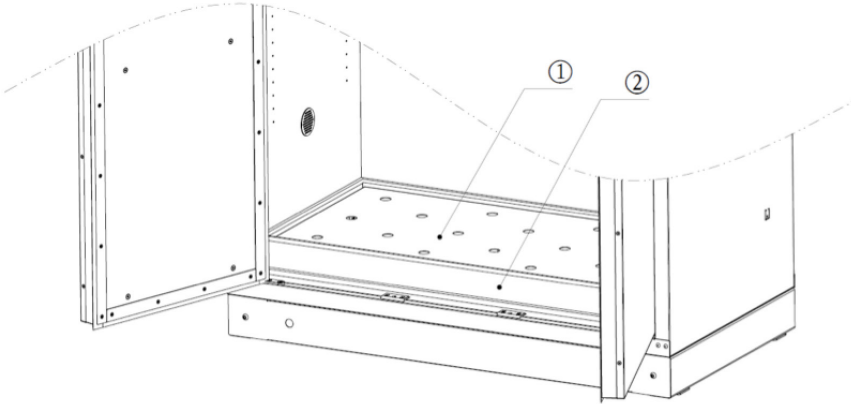
图 5 (picture 5)

i Attention

1. The cabinet door must be kept closed when not in use
2. Do not put your hands under the cabinet when lifting or moving the cabinet
3. The cabinet body must be placed horizontally to ensure that the cabinet door can be opened and closed normally
4. In order to ensure the air flow in the cabinet, please connect to a suitable ventilation device
5. Please place the cabinet for indoor use, not outdoors

C Trademark & Copyright

SYSBEL[®] is a registered trademark of SYSBEL, which is protected by law.

According to copyright law, it is forbidden to copy the whole or particle content of this manual without the written consent of SYSBEL. We have tried our best to ensure that the information in this manual is correct, and if the information is updated, we will not inform you.

You can check our official website at www.sysbel.com.cn any time.

☎ 86-4008-94-8808
🏠 www.sysbel.com

printed in shanghai 2022
Code No. SYS220719

SYSBEL®为SYSBEL 注册商标，受法律保护
SYSBEL® is the registered trademark of SYSBEL